


# CryoHolder



Instructions for use  
Technical sheet  
Logistic specifications  
Manufacturer information

# Instructions for use

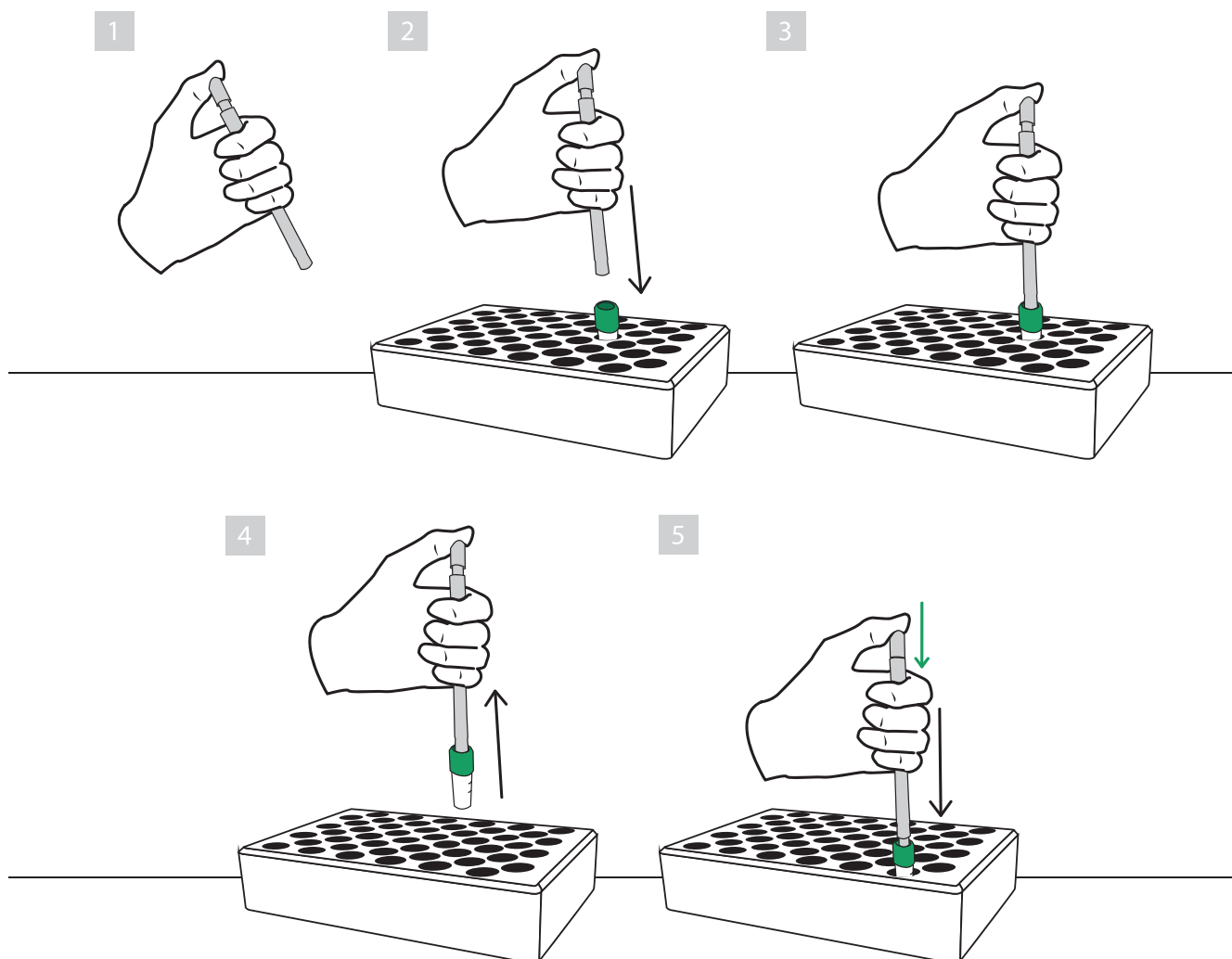

Prepare your CH for use.

1 Hold it in your hand with your thumb on the button at the top.

2 Choose the cryovial you want to move and point the CH at it.

3 Push the tip of the CH into the hole on the cap of the cryovial.

4 Now use the CH to move the vial to the new location.

5 Push the button with your thumb to release the vial.

Clean your CH after every use and store it properly.

Cryo  
Holder

**PharSol**

# Technical sheet



## CryoHolder

## Catalogue nr.: PS-001-CH

Product name:	CryoHolder
Dimensions :[mm]	141x10 (HxW)
Weight [g]:	200,00
Materials:	Casing and everything besides the spring - Aluminium 6026  Compression spring - stainless steel spring 1.43410 with standard EN 10270-3
Country of origin:	Slovenia, EU
Status of Intellectual property:	Patent Pending

### Description:

The Cryoholder is a tool for a safe, fast and easy transfer of frozen cryovials from one point to another. It is completely sterilisable in the autoclave and with disinfectant.

The tool is shaped like a pen and works similarly to a pen, with a button on the top and an inner piston which extends out of the tip of the tool.

### Main advantages:

- 100% safe transfer of cryovial
- Sterilisable
- Indestructible
- Easy-to-use
- Faster method of transfer

## Logistic specifications

M.O.Q.:	5 pcs
Products per:	
inner packaging:	28 pcs
pallet:	3584 pcs
Weight per:	
inner packaging:	2 kg
pallet:	256 kg + 25 kg
Customs Tariff Nr.:	HS Code 76169990
Delivery time:	5-30 days (depending on the quantity)

## Manufacturer information



+386 41 945 959  
info@pharsol.com  
www.pharsol.com

**Pharsol, d.o.o., Vreclova ulica 2, 2311 Hoče, Slovenia**  
VAT ID: SI25601580 | REGISTER NUM: 7196903000  
IBAN:SI56 1010 0005 5667 713 (INTESA SANPAOLO)

Cryo  
Holder

**PharSol**